

Closures - Cap Swabbing:



1 cap per 1 swab / 4 swabs needed

1. Remove swab from tube and swab the internal area of the cap.
2. Place swab back into the tube and return to the lab.




Closures - Cap submission:

1 cap per 1 swab / 4 caps needed

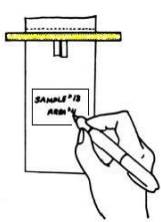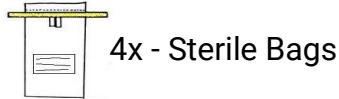


Complete label info.

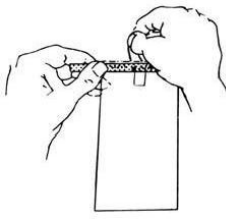


Remove top of bag at perforation.




Pull white tabs apart to open bag.

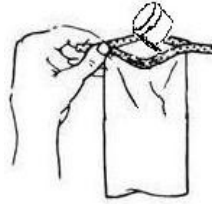

Insert 1 cap into bag. Do not touch inside of bag.

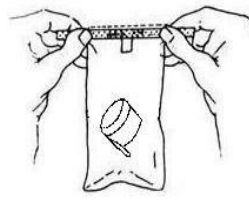


Roll over top of bag 3-4 times.

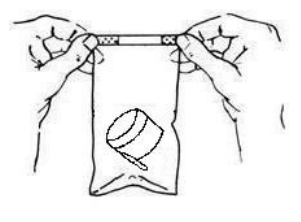

Fold in yellow tabs to lock bag

Containers

1 vial per container / 4 containers needed

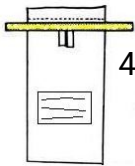


- Break Tamper Seal on bottle
- Twist off Lid.
- Do not touch contents inside surface of vial or lid.



Carefully empty contents of vial into container.



Shake and swirl to contact broth with entire inside surface area of container.

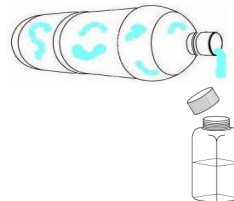

Carefully return liquid to original vial and close lid securely.



Insert each vial into its own sterile bag and close in the same manner as with the caps.

Shipping Instructions:

- Hold Time: 30 Hours
- Ship samples to: **Edge Analytical, 805 W. Orchard Drive, Suite 4, Bellingham, WA 98225** or drop off at lab.
- Make sure paperwork is properly filled out and **placed in a plastic bag** to protect it from being damaged.
- Ship "Next Day Air" / **Do Not ship samples on Thursday or Friday**

Place containers in cooler(s) with enough ice to keep samples at 4° Celsius +/- 2° until arrival at Edge.

- To prevent leaking, double bag the cubed ice.